

TruFidelis Plus DNA Polymerase

FB02005010 100 Reactions

FB02005050 500 Reactions

Store at -25°C to -15°C

TruFidelis Plus DNA Polymerase is a high-fidelity, proofreading DNA polymerase combining a novel enzyme with a processivity-enhancing domain. It features a hot start mechanism, allowing room temperature reaction setup and storage of pre-assembled PCR reactions. The polymerase is ideal for applications requiring high accuracy, such as cloning, sequencing, and site-directed mutagenesis.

- > 100X fidelity compared to Taq polymerase.
- 5'→3' DNA polymerase activity.
- 3'→5' exonuclease activity.
- Generates blunt-end amplification products.
- Amplifies up to 10 kb from genomic DNA and up to 20 kb from low complexity DNA.

Kit Contents

Reagents	100 Reactions	500 Reactions	Description
TruFidelis Plus DNA Polymerase	50 µL	250 µL	-
5X TruFidelis Plus Buffer	1.25 mL	5 x 1.25 mL	Provides 1.7 mM MgCl ₂ at 1X concentration
5X TruFidelis Plus GC Enhancer	1.25 mL	4 x 1.25 mL	Recommended for targets with >65% GC Sequence

PCR Reaction Setup:

For multiple reactions, prepare a master mix of components common to all reactions to minimize pipetting error, then dispense appropriate volumes into individual PCR tube before adding template DNA.

Component	20 µL Reaction	50 µL Reaction	Final Concentration
Water, nuclease-free	to 20 µL	to 50 µL	-
5X TruFidelis Plus Buffer	4 µL	10 µL	1X
10 mM dNTP Mix	0.4 µL	1 µL	0.2 mM each
Forward Primer	varies	varies	0.5 µM*
Reverse Primer	varies	varies	0.5 µM*
Template DNA	varies	varies	5-100 ng (gDNA) or 0.01-10 ng (plasmid DNA) per 50 µL reaction
TruFidelis Plus DNA Polymerase	0.2 µL	0.5 µL	-
5X TruFidelis Plus GC Enhancer (optional)	4 µL	10 µL	1X

*Reduce the final primer concentration to 0.2 µM when amplifying >5 kb targets from high complexity DNA and multiplex reactions.

Thermal Cycling Conditions on Thermal Cycler:

Step		2-step protocol (for primers >30 nt in length)		3-step protocol	
		Temp.	Time	Temp.	Time
Initial denaturation		98°C	30 sec	98°C	30 sec
25-35 PCR cycles	Denature	98°C	5-10 sec	98°C	5-10 sec
	Anneal	72°C	15-30 sec/kb	varies	10 sec
	Extend			72°C	15-30 sec/kb
Final extension		72°C	5 min	72°C	5 min
		4°C	hold	4°C	hold

Use your PCR product immediately in downstream applications, or store it at -20°C.